



**GAWANDE COLLEGE OF  
PHARMACY.**

Sakharkherda, Lavala Road, Tq.  
Sindkhed Raja  
Dist.Buldhana



Department: **Pharmaceutics**

Name of the instrument: **Ampoule Filling and Sealing Device**

S.O.P. No.: GCP/PH/035


Stock Register No.: PH-00106/79

Effective date: 27/2/14

Review date: 3/2/14

**Procedure:**

- Put plastic tube (right hand side bottom) in liquid vessel (or solution), is to fill in empty ampoule.
- Adjust the solution capacity by injection frame (working capacity 10ml to 50 ml).
- Keep the empty ampoule in front of center nozzle.
- Pull the injection inside & outside. The solution comes in ampoule.
- After filling the ampoule with solution, put the ampoule in hole of plate in the top of machine platform.
- Now rotate the ampoule with button provided in front of instrument and seal ampoule with gas burner.
- After use, clean injection valve and injection pump with distilled water.

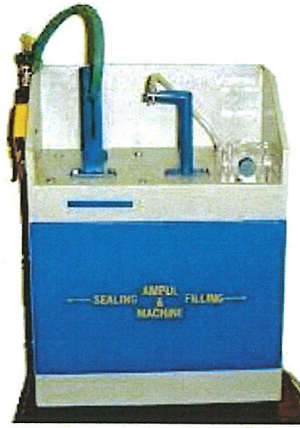

**Ampoule Filling and Sealing Device**

**Prepared by**

*S. Seelert*

Lab-In Charge  
*Miss S.D. Solat*

**Checked by:**

*Prof. A.G. Jadha*

**(Prof. A.G. Jadha)**  
HO.D

**Approved by**

*Dr. J.B. Sanap*

**(Dr. J.B. Sanap)**  
Principal