



**GAWANDE COLLEGE OF
PHARMACY.**

Sakharkherda, Lavala Road, Tq.

Sindkhed Raja

Dist.Buldhana



Department: Pharmaceutics

Name of the instrument: BULK DENSITY APPARATUS

S.O.P. No.: GCP/PH/038

**Stock Register No.: PH-
002167/01**

Effective date: 28/2/14

Review date: 30/2/14

Procedure:

- Weigh 20 g of product (to fill half of the cylinder) and transfer carefully in to the cylinder.
- Switch on the mains of the equipments.
- Put on the alarm switch.
- Put on the main of motors.
- After 50 strokes when a beeping sound is heard, put off the mains of the motor. Calculate the bulk density of the blend according to the formula.



BULK DENSITY APPARATUS

Prepared by

S. Sulek

Lab-In Charge

M.P.S.S. S.O.D. Sulek

Checked by:

A.G. Jadhav

(Prof. A.G.Jadhav)

H.O.D

Approved by

J.B. Sanap

(Dr..J.B.Sanap)

Principal