




**GAWANDE COLLEGE OF  
PHARMACY.**

Sakharkherda, Lavala Road, Tq.

Sindkhed Raja

Dist. Buldhana

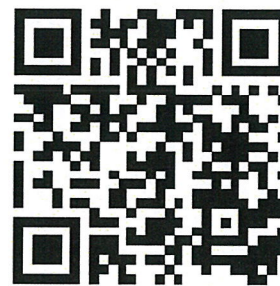

Department: **Pharmacology**

Name of the instrument: **CENTRIFUGE MACHINE**

S.O.P. No.: GCP/ PC/002

Stock Register No.: PH-002/02

Effective date: : 27/02/2014

Review date: 2/02/2014



**Procedure:**

1. Ensure that the instrument is clean and free from dust.
2. Ensure that all the knobs are in normal position.
3. Open the upper glass lid by releasing the lock and lifting it up.
4. Place the centrifuge tube in the compartment provided for it.
5. The compartment is design in such way to place 6 centrifuge tube at a time.
6. Switch 'ON' the mains.
7. Adjust the RPM speed by the knob provided on the front panel of the instrument.
8. A scale with deflection needle is provided on the front side panel of the instrument.
9. Set the required time by TIMER setting knob having range 0 to 60 minute.
10. Switch 'OFF' the mains when not required.

**Prepared by**

*J.M. Adhale*  
MISS. J.M. Adhale  
Lab-In Charge

**Checked by:**

*J.B. Sanap*  
**(Dr. J.B. Sanap)**  
H.O.D

**Approved by**

*J.B. Sanap*  
**(Dr. J.B. Sanap)**  
Principal