



**GAWANDE COLLEGE OF  
PHARMACY.**

Sakharkherda, Lavala Road, Tq.

Sindkhed Raja

Dist.Buldhana



Department: **Pharmacognosy**

Name of the instrument: **DISTILATION UNIT**

S.O.P. No.: GCP/PH/007

Stock Register No.: PG-  
0011/01

Effective date: 17/02/2015

Review date: 21/02/2015

**Procedure:**

1. Start the DM water line and check the conductivity of the of the DM water to be fed to the distillation still.
2. Fill the DM water surge tank completely with DM water.
3. Check out both the steam valve on the top of each column are open, if not open them.
4. Open the main steam inlet valve and let steam in to the column, close the drain valve.
5. Put 'ON' the main switch of the distillation steel panel and start the feed water pump.



**DISTILATION UNIT**

<p><b><u>Prepared by</u></b> <i>A. Ingle</i> <b>Mr. A.D. Ingle</b> Lab-In Charge</p>	<p><b><u>Checked by:</u></b> <i>RU</i> <b>Prof. R.R. Upma</b> H.O.D</p>	<p><b><u>Approved by</u></b> <i>J.B. Sanap</i> <b>(Dr..J.B.Sanap)</b> Principal</p>
--	---	---