



**GAWANDE COLLEGE OF
PHARMACY.**

Sakharkherda, Lavala Road, Tq.
Sindkhed Raja
Dist.Buldhana



Department: **Pharmaceutics**

Name of the instrument:HARDNESS TESTER

S.O.P. No.: GCP/PH/011

Stock Register No.:PH-0020/02

Effective date:17/02/15

Review date:19/02/15

Procedure:


1. Ensure that the instrument is clean and free from any residue of dust or tablet particle.
2. Hold the tablet between the jaw and nozzle in edgewise position.
3. Adjust the scale (by sliding), so that the zero on the scale coincides with the pointer.
4. Turn the screw knob slowly till the tablet breaks. The pressure indicated on the dial is in Kilogram per sq.cm.




Prepared by

Miss B.D. Solat
S. Seelot
Lab-In Charge

Checked by:


(Prof. A.G.Jadhao)
H.O.D

Approved by


(Dr..J.B.Sanap)
Principal