



**GAWANDE COLLEGE OF
PHARMACY.**

Sakharkherda, Lavala Road, Tq.

Sindkhed Raja

Dist.Buldhana



Department: **Pharmaceutical chemistry**

Name of the instrument: **Heating mantal**

S.O.P. No.: GCP/PH/028

Stock Register No.: PC 00323/04

Effective date: 17/2/15

Review date: 19/2/15

Procedure:

1. Place the flask containing sample to be heated on the mantle.
2. Switch on the heating mantle by connecting the power supply.
3. Adjust the desire temperature by using the temperature knob.
4. Switch OFF the power supply after desired time period.
5. Carefully remove the flask.

Precaution:

1. Avoid the contact with hot surface.
2. Avoid the over-heating.
3. Keep the surface clean.



Heating mantle

<p><u>Prepared by</u></p> <p><i>(Signature)</i></p> <p>Lab-In Charge <i>Miss. A. G. Kakde</i></p>	<p><u>Checked by:</u></p> <p><i>(Signature)</i></p> <p><u>(Prof. T.R. Sheikh)</u></p> <p>H.O.D</p>	<p><u>Approved by</u></p> <p><i>(Signature)</i></p> <p>(Dr. J.B. Sanap)</p> <p>Principal</p>
--	--	--