




**GAWANDE COLLEGE OF
PHARMACY.**

Sakharkherda, Lavala Road, Tq.

Sindkhed Raja

Dist. Buldhana

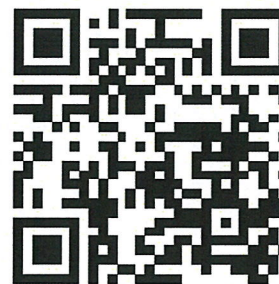

Department: **Pharmaceutical chemistry**

Name of the instrument: **Water Bath**

S.O.P. No.: GCP/PH/025

Stock Register No.: PC-0026/08

Effective date: 28/02/14

Review date: 2/03/14

/

Procedure:

1. Ensure the water level.
2. Connect the plug in the socket.
3. Switch ON the mains.
4. Adjust the required temperature using the setting knob.
5. Place the sample containing flask on the water bath station, for time specified in the procedure.
6. Switch OFF the mains.
7. Carefully remove the flask.


Precaution:

1. Avoid the contact with hot surface.
2. Avoid the over-heating of the sample.
3. Keep the surface clean.




Water Bath


Prepared by

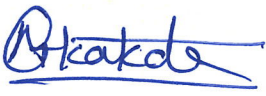
Miss. A. G. Kakade
Lab-In Charge

Checked by:


(Prof. T.R. Sheikh)
H.O.D

Approved by


(Dr. J.B. Sanap)
Principal